

## Tablet Triturate Mold

Hard-Coated (100 & 300 Cavities)

Two-piece molds designed for compounding tablet triturates and troches. Each mold consists of a base plate with alignment pegs and a top plate with corresponding holes. To make tablets, place the top plate on an Ointment Slab, Glass (SKU: E0595-00) or a glass pill tile, then fill the holes evenly with tablet paste using a hard rubber spatula to compress the mixture. When the paste has hardened, slide the mold off the edge of the slab, align the top plate alignment holes (left and right sides) over the base plate alignment pegs (left and right sides). Press down to punch out the tablets from the top plate. Allow the tablets to dry and harden.

Made of durable, corrosion and wear-resistant hard-coated aluminum, the hard coat is far superior to other "anodized" molds on the market in terms of scratch resistance, durability, and general use surface wear. Dishwasher safe.

### Tablet Triturate Mold

(100 Cavities)

SKU	Size	Qty
E0629-00	60 mg	Each
E0630-00	100 mg	Each
E0631-00	200 mg	Each

All are 6" x 4.5" / 15.2 cm x 11.4 cm.

### Tablet Triturate Mold

(300 Cavities)

SKU	Size	Qty
E0637-00	60 mg	Each
E0638-00	100 mg	Each
E0639-00	200 mg	Each

All are 6" x 11.75" / 15.2 cm x 29.8 cm.

	60 mg	100 mg	200 mg
Hole Size (Ø x D):	0.18" x 0.13" / 4.6 mm x 3.2 mm	0.24" x 0.13" / 6.0 mm x 3.2 mm	0.24" x 0.25" / 6.0 mm x 6.4 mm
Hole Volume:	0.053 mL	0.090 mL	0.181 mL

#### Recommended Cleaning Method:

Clean by hand with **Relianox™ Critical Cleaning Liquid Detergent** (SKU: P3023).

Hard-coated aluminum may be cleaned in an automatic dishwasher using mild detergents. Not recommended to use heavy acidic or caustic cleaning solutions. Do not clean with abrasive cloth, gritty materials, or steel wool. Use only plastic brushes or nylon/non-abrasive pads. Isopropyl Alcohol Cleaning & Drying: May be cleaned with isopropyl alcohol. Wipe with a cloth dipped in isopropyl alcohol, then immediately wipe dry with a clean cloth. May be dried in a hot air oven or using a portable hair dryer.

Autoclave safe. May be slight color variations from lot to lot.

#### See Also:

### TritTab™ SF/LF

Tablet Triturate Base

SKU	Size	NDC/Reference No.
B0215-40	100 g	72467-0215-40
B0215-50	500g	72467-0215-50
B0215-60	2.5 kg	72467-0215-60
B0215-65	5 kg	72467-0215-65

Specifically engineered for molded tablet triturates with the ability to use alcohol alone as the wetting solution and quickly disintegrate in the presence of moisture.

### TritTab™ Peppermint Ice

Natural Liquid Flavor Concentrate

SKU	Size	NDC/Reference No.
F0302-80	500 mL	72467-0302-80
F0302-85	3.8 L	72467-0302-85

Formulated specifically to be used alone as the wetting agent when compounding tablet triturates with **TritTab™ SF/LF Tablet Triturate Base** (SKU: B0215). Formulated with natural peppermint extracts. Water-soluble. 24-month BUD from date of manufacture.

This flavor can also be used as an effective water soluble liquid flavor concentrate. Suggested starting use level 0.1-0.5%.